

HOW TO CONNECT A GLOVE TO AN ISOLATOR SLEEVE - INSTRUCTION SHEET

Please use the instructions below as a guide to attaching a glove (e.g BFAP) to an isolator sleeve (e.g. GSL33NITPP26) with a BBCO-100 connector.



Prepare the isolator sleeve and connector ready for assembly, with the sleeve inside out



Take the sleeve and slide the connector over the bottom of it



Now, pull the bottom of the sleeve over the connector and tuck under the connector so that the sleeve is secured in place



Take the glove and feed it through the cuff connector, fingers first



Pull the cuff of the glove over the connector



Continue as in photo #5, ensuring to completely cover the sleeve cuff connector with the cuff of the glove



Take the silicone rubber O-ring and fix into groove of connector over the glove



Turn the assembled sleeve and glove the correct way round



Finally, place hand into fully assembled sleeve/glove system



